



Zymolyase[®] (from *Arthrobacter luteus*)

Zymolyase[®], produced by a submerged culture of *Arthrobacter luteus*⁽¹⁾, has strong lytic activity against living yeast cell walls^{(2),(3)} to produce protoplast or spheroplast of various strains of yeast cells. An essential enzyme for the lytic activity of Zymolyase[®] is β -1,3-glucan laminaripentaohydrolase. It hydrolyzes linear glucose polymers with β -1,3-linkages and releases specifically laminaripentaose as the main and minimum product unit^{(4), (5), (10), (11)}.

There are two preparations of Zymolyase[®], Zymolyase[®]-20T and Zymolyase[®]-100T, having lytic activity of 20,000 units/g and 100,000 units/g respectively. Zymolyase[®]-20T is ammonium sulfate precipitate while Zymolyase[®]-100T is a further purified preparation by affinity chromatography⁽⁹⁾. Lytic activity varies depending on yeast strain, growth stage of yeast, or cultural conditions⁽⁶⁻⁸⁾. Further informations related to Zymolyase[®] can be obtained in the reference section below⁽¹²⁻¹⁶⁾.



Specifications

Product Name		Zymolyase [®] -20T	Zymolyase [®] -100T
Form		Lyophilized Powder	
Purification		Ammonium Sulfate Precipitation	Affinity Chromatography
Activity		20,000 units/g	100,000 units/g
Essential enzyme		β -1,3-glucan laminaripentaohydrolase	
Other activities contained ^(*)	β -1,3-glucanase	approx. 1.5×10^6 units/g	approx. 1.0×10^7 units/g
	protease	approx. 1.0×10^4 units/g	approx. 1.7×10^4 units/g
	mannanase	approx. 1.0×10^6 units/g	approx. 6.0×10^4 units/g
Contaminants	Amylase, Xylanase, Phosphatase	Trace amount	Non detectable
Optimum pH and Temp.		pH7.5, 35°C (for lysis of viable yeast cells) pH6.5, 45°C (for hydrolysis of yeast glucan)	
Stability	2°C	No loss of activity was found after storage for 1 year	
Heat stability	30°C	70% of the lytic activity is lost after storage for 3 months	90% of the lytic activity is lost after storage for 3 months
	60°C	Lytic activity is lost on incubation for 5 minutes	
Specificity (Lytic Spectrum)		<i>Ashbya</i> , <i>Candida</i> , <i>Debaryomyces</i> , <i>Eremothecium</i> , <i>Endomyces</i> , <i>Hansenula</i> , <i>Hanseniaspora</i> , <i>Kloeckera</i> , <i>Kluyveromyces</i> , <i>Lipomyces</i> , <i>Metschnikowia</i> , <i>Pichia</i> , <i>Pullularia</i> , <i>Torulopsis</i> , <i>Saccharomyces</i> , <i>Saccharomycopsis</i> , <i>Saccharomycodes</i> , <i>Schwanniomyces</i> , etc.	

(*1) See reference, Kitamura, K., Kaneko, T., Yamamoto, Y., *J. Gen. Appl. Microbiol.*, **18**, 57 (1972) as to the definition of each enzyme units.


Unit Definition

One unit of lytic activity is defined as that amount which indicates 30% of decrease in absorbance at 800 nm (A_{800}) of the reaction mixture under the following condition.

[Reaction Mixture]	
Enzyme solution	: 1 ml (0.05-0.1 mg/ml for Zymolyase [®] -20T) (0.012-0.024 mg/ml for Zymolyase [®] -100T)
Brewer's yeast cell suspension	: 3 ml (2 mg/ml)
1/15M Phosphate buffer	: 5 ml (pH7.5)
Distilled water	: 1 ml

After incubation for 2 hours at 25°C with gentle shaking, A_{800} of the mixture is determined. When 60% of A_{800} decrease, equivalent to 2 units, is observed in the reaction system, the brewer's yeast cells are completely lysed, namely 1 unit of Zymolyase[®] lyses 3 mg dry weight of brewer's yeast.

Electron microscopical photo of yeast cell

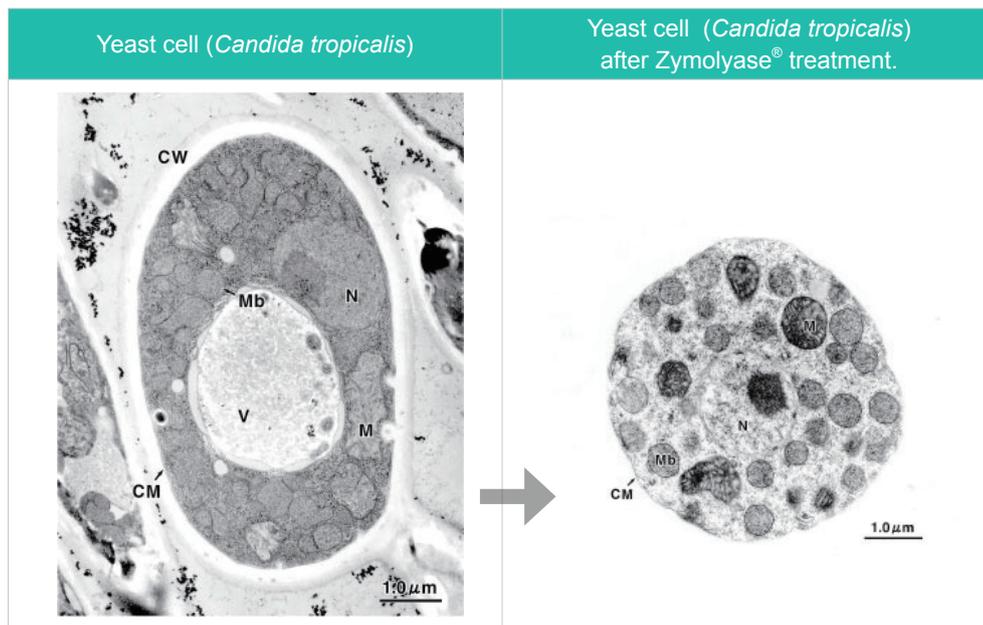

CW: Cell Wall
 CM: Cell Membrane
 M: Mitochondria
 Mb: Microbody
 N: Nucleus
 V: Vacuole

(Data courtesy of Masako Osumi, Emeritus Professor at Nippon Women's University)

Precaution on use

1. Avoid using nitrocellulose filters and use of material other than nitrocellulose when sterilizing. Zymolyase® may be adsorbed on nitrocellulose membranes.
2. Zymolyase® -100T may not be completely dissolved in buffers. Use Zymolyase® as suspension.
3. When sterilized, Zymolyase® is used in a concentration higher than 0.05%, prepare 2% Zymolyase® solution in buffers containing 5% glucose, filter the suspension and dilute the solution with the appropriate buffer.

Reference

1. Kaneko, T., Kitamura, K and Yamamoto, Y.: *J. Gen. Appl. Microbiol.*, **15**, 317 (1969)
2. Kitamura, K., Kaneko, T. and Yamamoto, Y.: *Arch. Biochem. Biophys.*, **145**, 402 (1971)
3. Kitamura, K., Kaneko, T. and Yamamoto, Y.: *J. Hen. Appl. Microbiol.*, **18**, 57 (1972)
4. Kitamura, K. and Yamamoto, Y.: *Arch. Biochem. Biophys.*, **153**, 403 (1972)
5. Kaneko, T., Kitamura, K. and Yamamoto, Y.: *Agric. Biol. Chem.*, **37**, 2295 (1973)
6. Kitamura, K., Kaneko, T. and Yamamoto, Y.: *J. Gen Appl. Microbiol.*, **20**, 323 (1974)
7. Kitamura, K. and Yamamoto, Y.: *Agric. Biol. Chem.*, **45**, 1761 (1981)
8. Katamura, K. and Tanabe, K.: *Agric. Biol. Chem.*, **46**, 553 (1982)
9. Katamura, K.: *J. Ferment. Technol.*, **60**, 257 (1982)
10. Kitamura, K.: *Agric. Biol. Chem.*, **46**, 963 (1982)
11. Kitamura, K.: *Agric. Biol. Chem.*, **46**, 2093 (1982)
12. Calza R. E., Schroeder A. L.: *J. Gen. Microbiol.*, **129**, 413 (1983)
13. Iizuka Masaru, Torii Yasuhiko, Yamamoto Takehiko: *Agric. Biol. Chem.*, **47** (12), 2267 (1983)
14. Shibata Nobuyuki, Kobayashi Hidemitsu, tojo Menehiro, Suzuki Shigeo: *Arch. Biochem. Biophys.*, **251**(2), 697 (1986)
15. Iijima Y., Yanagi S. O.: *Agric. Biol. Chem.*, **50** (7), 1855 (1986)
16. Herrero Enrique, Sanz Pascual, Sentandreu Rafael: *J. Gen. Microbiol.*, **133** (10), 2895 (1987)

Ordering Information

Product Name	Storage	Product No.	PKG Size
Zymolyase®-20T	R	07663-91	1 g
Zymolyase®-100T	R	07665-55	500 mg

Zymolyase® is a registered trademark of Kirin Holdings Company, Limited.

For research use only, not intended for diagnostic or drug use.

NACALAI TESQUE, INC.

Nijo Karasuma, Nakagyo-ku, Kyoto 604-0855 JAPAN

TEL : +81-(0)75-251-1730

FAX : +81-(0)75-251-1763

Website : www.nacalai.com

E-mail : info.intl@nacalai.com